

PRECAUTIONS : When using the long tip or extra-long tip

The liquid is affected by pressure, dependent on the size/length of the tip. Therefore when you use a certain pipette in combination with a long tip, it is necessary to adjust the scale on the pipette tip's height (=height of sample) in accordance to the level to which liquid has been drawn in.

If an 801* tip (250 μ L) or an 805** tip (1200 μ L) is used, it is necessary to make adjustments.

Note: The shape/size of the following tips is identical.

***Tip # 801: Tip # 200XS series**

****Tip # 805: Tip # 1200S series**